

Biological Microscope



CX3 Series

Comfortable, High-Throughput Routine Microscopy







Comfortable for Long Periods of Routine Microscopy

The CX33 and CX43 microscopes enable users to remain comfortable during routine microscopy. The microscope frame well fit the hands and the location of the control knobs maximize ergonomics to improve work efficiency. Users can quickly set a specimen with one hand, while adjusting the focus and operating the stage with the other hand with minimal movement. Both microscopes also feature a camera port for digital imaging.





Maintain Preferred Observation Conditions with Minimal Adjustments

Uniform Illumination with Consistent Color Temperature

The color temperature of the CX LED illumination produces daylight conditions, so specimens can be viewed with their natural colors. The color temperature is consistent at any brightness, so users don't have to spend time making adjustments when they change brightness. The LEDs have a long 60,000-hour lifetime in the design value, helping reduce cost, and the brightness level remains stable throughout the LED's life.



Excellent Optical Performance for Flat Images

The microscope employs Plan Achromat objectives, which provide clear images with high image flatness over a wide field of view. This helps users view specimens clearly and evenly during routine microscope observations.



Select and Set Your Contrast Level

Users can preserve their favorite contrast by locking the aperture diaphragm. It stays fixed at the optimally chosen position if it is accidentally touched while changing slides.



Change Magnification without Adjusting the Condenser

Users can change the magnification from 4X to 100X without moving the top lens on the condenser. 2X magnification is also available by simply setting the objective and the condenser turret to 2X position.



Simple Fluorescence Observation

Fluorescence observation is simple and easy. Plug the compact fluorescent illuminator into the microscope frame for fluorescence observation. Its LED light source is pre-centered, and the transmitted illumination is shuttered by simply setting the condenser turret to the FL position. This reduces background noise in the fluorescence image from incidental light coming from the top lens of the condenser.



Remain Comfortable during Extended Usage

Single-Handed Sample Placement

A specimen can be quickly slid in and out using one hand. The specimen holder opens a little and firmly retains the specimen during operation. The versatile holder accommodates a variety of slide types, including a hemocytometer.



Use Up to Five Objectives

For added flexibility, up to five objectives can be supported by the revolving nosepiece. In addition to general objectives, users can select a 2X objective for wide area observation or objectives for phase contrast. These objectives with long working distances help keep specimens from getting damaged.



Ergonomically-Positioned Focus Knob

The low-positioned focusing knob enables users to make observations while keeping their hands and forearms rested on the desk, helping provide comfort. The focusing stopper prevents a specimen from accidentally hitting an objective when working under high magnification.

Smooth Magnification Change

The low-positioned revolving nosepiece enables users to quickly change magnifications with minimal arm movement between focusing, greatly improving work efficiency during prolonged use.





Ergonomic Stage and Eyepiece Position

The low-positioned stage is designed to enhance comfort and reduce fatigue. The stage surface can be widely seen from the eye point position, which enables users to smoothly set and check specimens on the stage. The stage knob can be controlled with just a light touch and can be adjusted at the same time as the focusing knob, since they are located close together.



Specimen Holders that Match Your Observation Style

Stage accessories improve efficiency when users need to observe a large number of specimens. With the specimen holder sheet, a specimen can be freely operated by a finger on the sheet and can be precisely adjusted using the stage knob. The double specimen holder can retain a large specimen or two specimens.



Simplified Fluorescence Observation

Fluorescence observation can be easily set up on the standard configuration while keeping the eye point the same as other observation methods. Simply plug the compact fluorescent illuminator into the back of the microscope frame.



Versatile Applications

The universal condenser offers a variety of observation methods and future upgradability. In combination with the five-position revolving nosepiece, multiple applications can be covered using the single microscope frame.

Brightfield



Leukocyte (minimum iris aperture)

Brightfield

Fluorescence



Urinary Cast (minimum iris aperture)



HeLa cells

Accessories

Phase Contrast

Simple polarizing intermediate attachment/CX3-KPA

Offers polarized observation of urate crystals and amyloid in combination with a polarizer and analyzer.

Eyepoint adjuster/ U-EPA2

Raise the eyepoint position by 30 mm for added comfort.

Arrow pointer/ U-APT

Insert an LED arrow into your image; great for digital imaging and presentations.

Dual observation attachment/U-DO3

Enables dual, simultaneous observation of a single specimen from the same direction with equal magnification and brightness for both operators. A pointer can be used to indicate specific sections of the specimen to simplify the training process and enhance discussion.













Renal Glomerulus

CX33 Microscope

For less demanding requirements using only brightfield and darkfield, the CX33 microscope is a great option. The low-positioned nosepiece and stage, focusing lock, specimen holder, and inward quadruple revolving nosepiece make the CX33 microscope is well-suited for everyday observations in one easy configuration.



CX43 Specifications

Optical System		UIS2 (universal infinity-corrected) optical system							
Illumination System		Built-in transmited illumination system Köhler illumination (fixed field diaphragm) LED power consumption 2.4 W (nominal value), precentered							
Focusing		 Stage height movement (coarse movement stroke: 15 mm) Stroke per rotation for coarse adjustment knob: 36.8 mm, Focusing stopper Torque adjustment for coarse adjustment knob Fine focus knob (minimum adjutsment gradations: 2.5 µm) 							
Revolving Nosepiece		Fixed quintuple nosepiece with inward tilt							
Stage		 Wire movement mechanical fixed stage, (W × D): 211 mm × 154 mm Traveling range (X × Y): 76 mm × 52 mm Single specimen holder (optional: double specimen holder, sheet holder) Specimen position scale Stage XY movement stopper 							
Observation Tube	Type (anti-fungal)	Binocular		Trinocular			Tilting binocular		
	Eyepiece (anti-fungal)	10X Field Number (FN): 20		10X Field Number (FN): 20			10X Field Number (FN): 18		
	Tube Inclination	30°		30°			30°-60°		
	Light Path Selector	None		None (eyepiece/camera port = 50/50 fixed)			None		
	Interpupillary Distance Adjusting Range	48-75 mm							
		 Universal condenser with 7 turret positions: BF (4-100X), 2X, DF, Ph1, Ph2, Ph3, FL Condenser turret lock pin (BF only) Built-in aperture iris diaphragm AS lock pin 							
Observation Methods		Brightfield, simple polarization, fluorescence, phase contrast, darkfield							
Objectives		Plan achromat (UIS2 2X NA 0.0 4X NA 0.1 10X NA 0.2 20X NA 0.4 40X NA 0.4 60X NA 0.5 100XOO NA 1.2 100XOI NA 1.2	2), anti-fungal D6 W.1 1 W.1 25 W.1 4 W.1 35 W.1 35 W.1 25 W.1 25 O.6 W.1	 D. 5.8 mm D. 18.5 mm D. 10.6 mm D. 1.2 mm D. 0.6 mm D. 0.2 mm D. 0.13 mm D. 0.13 mm 	10XPH 20XPH 40XPH 100XOPH	NA 0.25 NA 0.4 NA 0.65 NA 1.25	5 W.D. W.D. 5 W.D. 5 W.D.	10.6 mm 1.2 mm 0.6 mm 0.15 mm	
Fluorescence Light Source		Easily add an LED reflected fluorescence illuminator (peak excitation wavelength 470 nm: B excitation only), precentered							
Rated Voltage/Electric Current		AC 100-240 V 50/60 Hz 0.4 A							
CX33 Spec	ifications								
Optical System		Infinity optical system							
Illumination System		 Built-in transmited illumination system Köhler illumination (fixed field diaphragm) LED power consumption 2.4 W (nominal value), precentered 							

	LED power consumption 2.4 w (nominal value), precentered					
Focusing	Stage height movement (coarse movement stroke: 15 mm) Stroke per rotation for coarse adjustment knob: 36.8 mm, Focusing stopper Torque adjustment for coarse adjustment knob					
	 Fine focus knob (minimum adjutsment gradations: 2.5 µm) 					
Revolving Nosepiece	Fixed quadruple nosepiece with inward tilt					
Stage	 Wire movement mechanical fixed stage, (W × D): 211 mm × 154 mm Traveling range (X × Y): 76 mm × 52 mm Single specimen holder (optional: double specimen holder, sheet holder) Specimen position scale Stage XY movement stopper 					
Observation Tube	 · 30° inclined trinocular tube (anti-fungal) · Light path selector: eyepiece/camera port = 100/0 or 0/100 · Interpupillary distance adjusting range: 48-75 mm · Eyepoint adjustment: 375.0-427.9 mm 					
Eyepieces (anti-fungal)	· 10X Field Number (FN): 20 · 15X Field Number (FN): 16 (optional)					
Condenser	Abbe condenser NA 1.25 with oil immersion Built-in aperture iris diaphragm					
Observation Methods	Brightfield, darkfield					
Objectives	Plan achromat, anti-fungal 4X NA 0.1 W.D. 27.8 mm 10X NA 0.25 W.D. 8.0 mm 20X NA 0.4 W.D. 2.5 mm (optional) 40X NA 0.65 W.D. 0.6 mm 100X NA 1.25 W.D. 0.13 mm (optional)					
Rated Voltage/Electric Current	AC 100-240 V 50/60 Hz 0.4 A					

OLYMPUS CORPORATION is IS014001 certified.
 OLYMPUS CORPORATION is IS09001 certified.
 OLYMPUS CORPORATION is IS013485 certified.

All company and product names are registered trademarks and/or trademarks of their respective owners.
 Specifications and appearances are subject to change without any notice or obligation on the part of the manufacturer.

www.olympus-lifescience.com

OLYMPUS CORPORATION Shinjuku Monolith, 2-3-1 Nishi-Shinjuku, Shinjuku-ku, Tokyo 163-0914, Japan